Highlighting innovative design features and useful application information for


Thermo Scientific CO₂ Incubators.

smart

design and innovation CO₂ Incubation





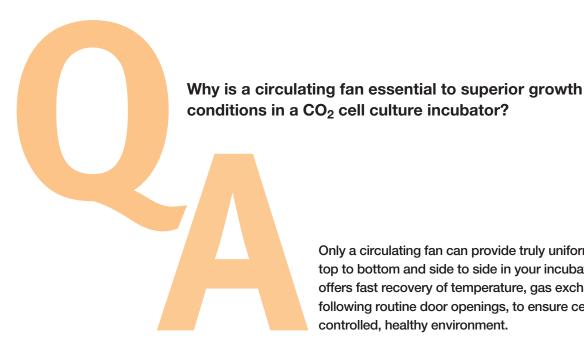

Only a circulating fan can provide truly uniform conditions from top to bottom and side to side in your incubator. A circulating fan offers fast recovery of temperature, gas exchange and humidity, following routine door openings, to ensure cells grow in a well controlled, healthy environment.

When any incubator door is opened, physics dictate that the conditions inside rush to equilibrate with the external atmosphere. Once the door is closed, the use of an integrated fan provides the fastest recovery to your desired conditions, thus minimizing stress to cells from loss of temperature, CO2 and humidity



Why Thermo Scientific CO₂ Incubators?

Active airflow circulation provides the most uniform growth conditions

Without a uniform environment throughout your CO_2 incubator, cells in different areas can experience varying conditions, leading to confusing results. Thermo Scientific fan-assisted incubator designs ensure homogeneous conditions throughout the entire chamber, so that all cells experience the same temperature, gas exchange and humidity, regardless of their location in the growth chamber. These standardized conditions form the basis of good, controlled science and meaningful, reproducible results.

Without a fan, stratification of critical incubator conditions can compromise research results

An incubator with no fan relies upon gravity convection, or slow moving thermal currents, for humidity transport and gas exchange. This design results in slow recovery rates such that frequent door openings can prevent your cultures from experiencing adequate exposure to your selected culturing conditions. Lack of uniformity within the chamber can create variation among your cell cultures with a negative impact upon your project goals. This is especially critical in an incubator with variable oxygen control, since oxygen concentration affects nearly every cellular process.


Incubators without fans do not benefit from the positive pressure that helps to prevent the entry of unwanted environmental microorganisms during door openings. The airflow systems in Thermo Scientific incubators are specifically engineered to provide a continuous, gentle flow pattern directed around the sides of the chamber to prevent disruption, evaporation and desiccation of culture medium.

Summary


Thermo Scientific CO₂ incubators with precisely designed fan assisted airflow systems provide more efficient circulation of critical temperature, gas and humidity conditions to provide healthy, uniform cells.

Comparative Recovery Rates of Thermo Scientific Fan-assisted CO₂ Incubators vs. Fanless Designs


Temperature Recovery Rates

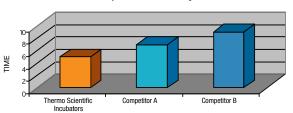

CO2 Recovery Rates

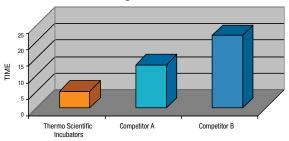

rH Recovery Rates

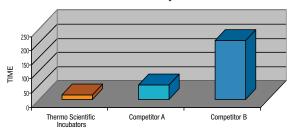



Note: Data was collected following a 30 second door opening, 5% CO₂, 22°C ambient. Results amay vary with operating conditions.

See how Thermo Scientific CO₂ incubators provide optimal cell growth. Learn more at www.thermoscientific.com/co2incubators

© 2015 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries.

Australia +61 39757 4300 Austria +43 1 801 40 0 Belgium +32 53 73 42 41 China +800 810 5118 or +400 650 5118 France +33 2 2803 2180

Germany national toll free 0800 1 536 376 Germany international +49 6184 90 6000 India toll free 1800 22 8374 India +91 22 6716 2200 Italy +32 02 95059 552 Japan +81 3 5826 1616 Netherlands +31 76 579 55 55 New Zealand +64 9 980 6700 Nordic/Baltic/CIS countries +358 9 329 10200 Russia +7 812 703 42 15 Spain/Portugal +34 93 223 09 18 Switzerland +41 44 454 12 22 UK/Ireland +44 870 609 9203 USA/Canada +1 866 984 3766

Other Asian countries +852 2885 4613 **Countries not listed** +49 6184 90 6000

